

photometrics®  
**CoolSNAP™**  
**MYO**


1940 x 1460 imaging array  
4.54 x 4.54 μm pixels

The CoolSNAP MYO is a high resolution, high sensitivity camera for moderate to low-light life science applications. This unique cooled CCD provides 4.54μm pixel pitch, 14-bit digitization at 20MHz, enabling high spatial resolution and an optimized frame rate for live cell imaging. Its 2.8 Megapixels and a high Quantum Efficiency enables sensitive imaging with the option for binning for a higher dynamic range as well as increased signal-to-noise performance – all while providing an ideal pixel pitch for microscopy.

**Primary applications**

- Fixed Cell Imaging**
- Immunofluorescence**
- Cell Trafficking**
- FRET, FRAP, FISH**
- Near-Infrared DIC**
- Calcium/Ion Imaging**

Features	Benefits
1940 x 1460 imaging array 4.54 x 4.54 μm pixels	High spatial resolution for imaging finer details
High Quantum Efficiency	~75% peak quantum efficiency delivers high sensitivity
20 MHz read out	High Speed readout to maximize temporal resolution
USB 2.0 Interface	Easy connectivity and setup
Binning	Increase frame rate and signal-to-noise performance
14-bit digitization	Quantify bright and dim signals in the same image
Thermoelectric cooling	Stabilized cooling produces a low dark current for long exposures
Fan Disable Option	Disable the fan for vibration-sensitive applications
C-mount	Easily attaches to microscopes, standard lenses, or optical equipment
Acquisition software	Captures, analyzes, and saves high-resolution images
PVCam® Driver	Support in a wide range of third party software packages Supported in Windows 7 64-bit/32-bit

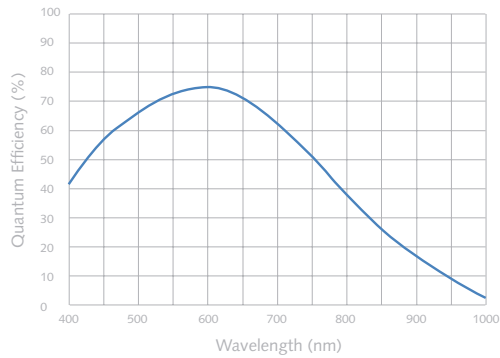


Binning	Region		
	1940 x 1460	970 x 730	646 x 486
1 x 1	6.3	11.8	16.8
2 x 2	11.6	20.8	28.4
3 x 3	16.2	27.8	36.5
4 x 4	20.1	33.2	42.6

(Frames per second)

Note: Frame rates are measured at 20 MHz with 0-millisecond exposure times.

Specifications	
CCD Sensor	Sony® ICX-674 Interline CCD
CCD Format	1940x1460 imaging array 4.54x4.54 μm pixels 8.8 x 6.6 mm imaging area (11mm diagonal, 2/3" format)
Linear Full-Well	12,000e-
Read Noise	5.5e- @ 20MHz 4.5e- @ 10MHz 3.5e- @ 1.25MHz
Nonlinearity	<1%
Digitization	20MHz, 10MHz, 1.25MHz
Cooling	0°C
Dark Current	0.005 e-/pixel/second @ 0°C
Operating Environment	0 to 30°C ambient, 0-80% relative humidity non-condensing
Triggering	Trigger First Mode Strobe Mode Bulb Mode
Power Requirements	5V DC, 4A Maximum

Note: Specifications are typical and subject to change.

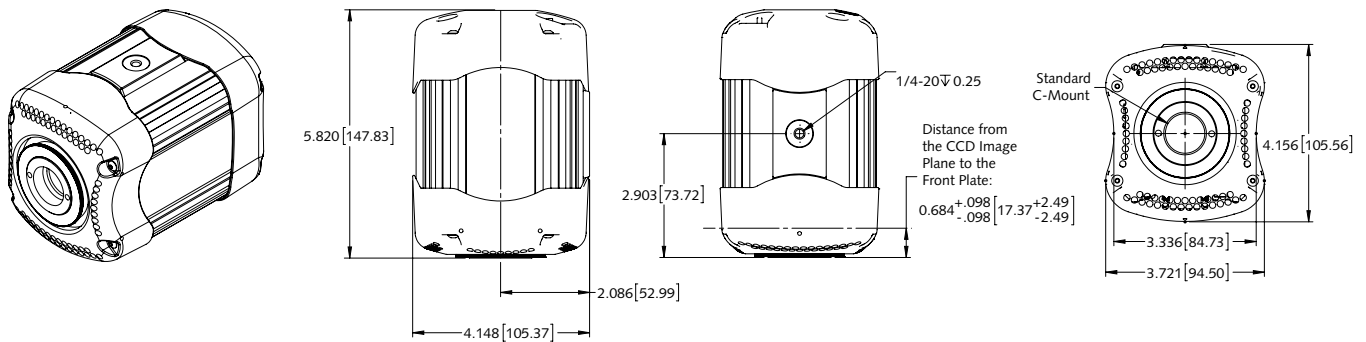

CoolSNAP is a trademark of Photometrics. Photometrics and PVCam are registered trademarks of Photometrics. Other brand and product names are the trademarks or registered trademarks of their respective owners and manufacturers.

USA 520.889.9933  
Asia Pacific +65.6841.2094

France +33.1.60.86.03.65  
Germany +49.89.660.779.3

Japan +81.3.5639.2731  
UK +44.1628.890858

**PHOTOMETRICS®**  
www.photometrics.com  
info@photometrics.com